

## VENUS ONE service pack 1

The service pack 1 CD contains the following main changes:

### Enhancements:

- ▶ Maintenance methods for the Nano pipettor.
- ▶ Maintenance methods for the 5ml channels.
- ▶ Optimized liquid following in mix mode independent from Asp/Disp volume.

### Fixes for the following issues:

- ▶ Optimized Report Mapping File using needles with mix settings (SCR 1385).
- ▶ Single Image Channels support (SCR 1386).
- ▶ Support for dynamically added labware (SCR1394).
- ▶ Position correction of CO-RE 384 Head using RocketTips on -X (SCR 1398).
- ▶ Method editor closes when ending Run Control with open workflow (SCR 1400).
- ▶ Support for sequences of attached system decks in step combo box (SCR 1402).
- ▶ Support for file selection dialogs under Windows VISTA (SCR1404).
- ▶ Implementation of the Software Change Requests 1405 and 1407.

### Installation:

The installation of this service pack cd can be performed over an existing VENUS one version. It is not necessary to uninstall the installed version.

Only MICROLAB STAR instruments with the latest firmware package 5.7S (including the master firmware 6.1S) will be supported by this service pack.

This is due to the integration of the enhanced mixing liquid following algorithm for the MICROLAB STAR pipetting components like pipetting channels or heads.

