

HAMILTON METHOD MANAGER v5 – Alvaro Cuevas 2010

**This new version is compatible with Vista and Windows 7, and has a simulation toggle button.

This program is created to help you to organize your Hamilton methods and speed up the launch process. The full application is designed to be easily adapted to new touch screen PCs.



Methods are packed in “groups of methods”. You can select the icon, name and color of your group to easily recognize them in the screen.

You can link to each method files, pictures, or any additional info you need.

New to Version 5: HamiltonPath text file

Please write the path to your HAMILTON folder in the HamiltonPath.txt file. This file has to be in the same folder that the MethodManager.exe and the Manager_DB.mdb.

Remember to end the path with a backslash “\”, and please don’t use “” at all.

E.g. c:\Program Files\Hamilton\


There are 2 view modes in the application, you can switch modes clicking button at the top right corner.

- **Run Mode:** Default. Run methods, view all the info and method attachments. At the bottom of this screen also appears “Quick buttons” to the Frequent Used Methods (method launched in HxRun Control from this app).



Just click one group (big squares) to see the related methods, and click it again to close it and back.


You can switch simulation mode ON or OFF with the button at the top right

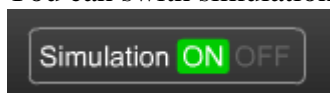



Clicking the “Run” button of each method, will open it in the Hamilton Run Control.
 Clicking “Layout” or “Anex” will open the linked file with the default system application.
 Clicking the “Summary” will pop up the short description of the method.

- **Edit Mode:** Add / Delete groups and methods.

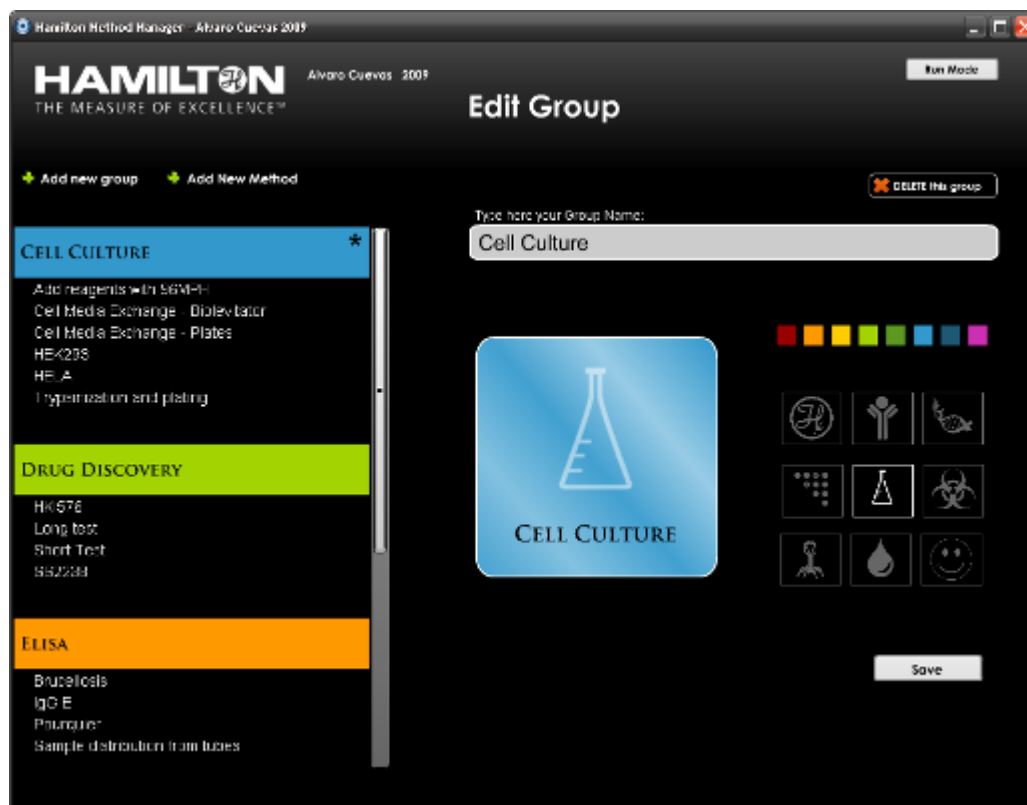


In the edit mode, click buttons **+ Add new group** **+ Add New Method** at the top left to add new items.

For editing existing items, just click on the left menu.

If you click a group (colored header title) the edit group screen will show (See previous screenshot). From here you can change color, icon, name of the selected group. Just remember to click “SAVE” button at bottom to save changes. You can delete the group clicking the button at the top of the Name input field.

If you click a method, the edit method screen will show. Here you can link the *.med, *.hsl or *.wfl file, layout pictures, videos, pdf, doc or any file you want to attach to the method profile.

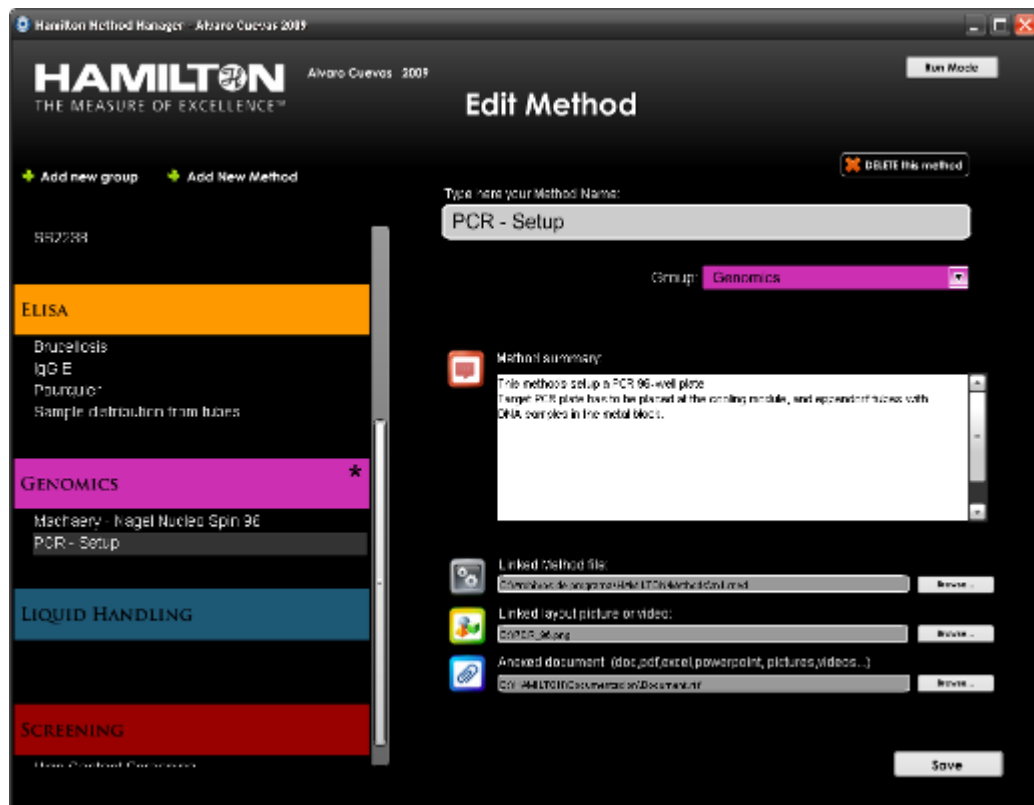

New is in this version:

- Recoded function that automatically resizes the group title as you type, to work on new computers.
- Restricted the text input to not allow quotes (") that were crashing the Access database of the Method Manager in some computers.

om

Run Mode

Edit Group

✗ DELETE this group

Type here your Group Name:

group of methods of my crazy customer



GROUP OF METHODS OF MY CRAZY CUSTOMER



















Group name should be automatically resized to fit in here as you type it above

Save