

## **MICROLAB® STAR VENUS two Software – Service Pack 1**

Date: 2011-05-16 (Build 4785)

**This service pack contains the following main changes:**

- **Enhancements:**
  - Maintenance methods for CO-RE 96 TADM head.
- **Fixes for the following issues:**
  - Using Custom Dialog Step within workflow (impacts only systems using VENUS Dynamic Scheduler).
  - Container z-height values for irregular rack definition.
  - Liquid class parameter within ML STAR Steps (use of arrays).
  - Sequence view optimization when using probe stamp for highlight.
  - Compatibility to existing drivers (e.g. HxFan, ELx405).
  - Use of the “Control Panel-Initialize” to eject tips after method abort.

**Please note that it is not possible to execute a “repair installation” on the basepack when the service pack is installed. Both the service pack and base package have to be removed and reinstalled if a “repair installation” is needed.**