



# VENUS 6.0.1

Release notes | 2023-06-09

---

## STAR and VANTAGE:

Function	Description / Changes / New features
<b>General</b>	<ul style="list-style-type: none"><li>• Support of Windows 10 64-bit and Windows 11 64-bit operating system</li></ul>
<b>Phoenix</b>	<ul style="list-style-type: none"><li>• Use of Phoenix 6.0.1.50042</li></ul>
<b>Interface</b>	<ul style="list-style-type: none"><li>• New application interface (single desktop icon)</li><li>• Dark and Light Theme options</li><li>• New - Home section with instrument status, recent runs and frequent shortcuts</li><li>• New - Shortcuts section</li><li>• New - Shortcut validation flag</li><li>• New - Run Control</li><li>• New - Trace Viewer</li><li>• New - Run History section</li><li>• New - 'Generate Diagnostics File' from run, with traces and method pkg</li><li>• New - Maintenance section</li><li>• New - Instrument Control section</li><li>• New - System Tools section</li><li>• New - Settings section with application settings and shortcuts manager</li><li>• New - Import/Export shortcuts and methods function</li><li>• New - Run History auto-cleanup option</li><li>• New - Support section for contact cards</li></ul>
<b>Remote Monitoring</b>	<ul style="list-style-type: none"><li>• New remote run monitoring over the web</li></ul>
<b>Editors</b>	Method Editor: <ul style="list-style-type: none"><li>• New - Add/remove multiple dependencies at once</li><li>• New - Add/remove multiple libraries at once</li><li>• New - Complex 'IF/ELSE' command with AND/OR</li><li>• New - Complex 'Assignment with calculations' command</li><li>• New - Support for emailing with SSL and TLS</li><li>• New - Simulation toggle under Tools menu</li></ul>

	<ul style="list-style-type: none"> <li>• New - Stop timer command</li> <li>• New - Stop individual timers at runtime</li> <li>• New - Timer display name</li> <li>• New - Sequence : Get remaining positions command</li> <li>• New - Validate method and all related files (except liquid classes)</li> <li>• New - Variables view context menu : Remove unused</li> <li>• New - Search with regular expressions, find previous, mark all</li> <li>• New - Ungroup steps</li> <li>• New - Group description (auto traced) and notes (not traced)</li> <li>• New - Trace timers start/end</li> <li>• New - Add multiple libraries and remove all libraries</li> <li>• New - Key shortcuts to group/ungroup</li> <li>• New - Allow editing a method while it's running (requires restart the run to apply changes)</li> </ul> <p>System Deck Editor:</p> <ul style="list-style-type: none"> <li>• Sequences list - highlight item on key press</li> </ul> <p>Labware Editor:</p> <ul style="list-style-type: none"> <li>• New - Support for Multiflex Carrier definition</li> <li>• New - Labware category manager</li> <li>• New - Labware filters</li> <li>• New - Automatic labware index generation</li> </ul> <p>Liquid Class Editor:</p> <ul style="list-style-type: none"> <li>• Support for MagPip liquid classes</li> </ul> <p>System Configuration Editor:</p> <ul style="list-style-type: none"> <li>• New - VANTAGE arms configurator</li> <li>• New - Email SMTP security protocol setting</li> <li>• Hamilton Authentication System option has been removed</li> </ul>
<p><b>Power Steps</b></p>	<p>The following Power Steps are available:</p> <ul style="list-style-type: none"> <li>• New - Load And Match</li> <li>• Transfer Samples</li> <li>• Add Reagent</li> <li>• Serial Dilution</li> <li>• Replicates</li> <li>• Hit Picking</li> </ul>
<p><b>Custom Dialog Step</b></p>	<ul style="list-style-type: none"> <li>• Included Template 'Hamilton'</li> <li>• Option to link external image</li> <li>• Default keep image aspect ratio</li> </ul>

	<ul style="list-style-type: none"> <li>• Define control tab index</li> <li>• Auto-close dialog after timeout</li> <li>• Multi-select controls</li> <li>• Align and distribute multiple controls</li> <li>• Option to set focus to specific control on startup</li> <li>• Startup dialog window location</li> <li>• Text input as password (mask input)</li> <li>• Text input automatic scrollbars</li> <li>• Use array values as in/out variable for checkboxes and radio buttons</li> </ul>
<b>Installer</b>	<ul style="list-style-type: none"> <li>• Activation of .NET 3.5 feature (internet access required) and check for .NET 4.6.2 availability</li> <li>• Microsoft® SQL Server 2019 Express Edition</li> <li>• Separate installers for STAR and VANTAGE</li> <li>• TADM installed by default (requires manual activation on STAR)</li> <li>• Scheduler installed by default</li> </ul>
<b>VANTAGE</b>	<ul style="list-style-type: none"> <li>• Support for dual arm configurations (no duplicated tools allowed)</li> <li>• New – MagPip Channels steps</li> <li>• New – 2T labware definitions for different cover types</li> <li>• New – Track Gripper startup maintenance routine</li> <li>• New – Track Gripper calibration alignment routine</li> <li>• New – 5mL daily maintenance routine</li> <li>• New – MagPip daily maintenance routine</li> <li>• Improved CAS (Collision Avoidance System)</li> <li>• Improved QCG error handling on abort</li> </ul>
<b>Bug Fixes</b>	<ul style="list-style-type: none"> <li>• Bug 34599: Status Light steps (Set and Progress) don't work when no Initialize step executed before</li> <li>• Bug 34808: Load Carrier - Barcode filename with parenthesis in the path cannot be used</li> <li>• Bug 36972: iSWAP does not re-open gripper when looking for the next plate on Get First Plate on Stack</li> <li>• Bug 17937: HxRun crash (exception code 0xc015000f) when calling SeqGetPositionCountForCurrLabware with no positions available</li> <li>• Bug 33304: Power Steps - Unnecessary tip pick up after processing all non-controlling seq positions</li> <li>• Bug 38453: Power Steps - Hit Picking - only first 8 volumes in the worklist are used</li> <li>• Bug 29443: Power Steps - Serial Dilution - Wrong volume in 2nd dilution if 'Pipette Samples' is selected</li> <li>• Bug 32077: Reading from *.xlsx file with no header skips first row</li> </ul>